tasco®

DIGITAL MICROSCOPE

INSTRUCTION MANUAL

MANUEL D'INSTRUCTIONS
MANUAL DE INSTRUCCIONES
BEDIENUNGSANLEITUNG
MANUALE DI ISTRUZIONI
MANUAL DE INSTRUÇÕES

Model#: 780200T

Lit#: 93-0599 / 12-10

Pollen	Apple	Hydrilla Leaf	Broad Bean Leaf	Onion Bulb Skin
Pollen	Pomme	Feuille d'Hydrilla	Feuille de fève	Pelure d'Oignon
Polen	Manzana	Hoja de Hydrilla	Hoja de haba	Piel de cebolla
Blütenstaub	Apfel	Blatt der Grundnessel	Blatt der Saubohne	Zwiebelhaut
Polline	Mela	Foglia di hydrilla	Foglia grande fagiolo	Pelle bulbo di cipolla
Pólen	Maçã	Folha da Hydrilla	Folha de Favas	Epiderme do Bulbo da Cebola

Sea Salt	Brine Shrimp Eggs	Gum Media	Eosin Dye
Sel de mer	Oeufs d'artemia	Milieu de gomme	Colorant éosine
Sal marina	Huevos de gamba en salmuera	Medio gomoso	Tinte de eosina
Meersalz	Salzwasser-Garneleneierl	Gum-Media	Eosin-Farbe
Sale marino	Uova di gambero in salamoia	Elemento in gomma	Colore eosina
Sal Marinho	Ovos de Artémia	Meio Resinoso (Bálsano do Canadá)	Corante Eosina

TABLE OF CONTENTS

Content	Page
Safety Information	4
Technical Data	4
System Requirements	5
Included Items	5
Parts Guide	6
Assembly	7
Software Installation	7
Selecting the Lighting	8
Setting the Magnification	8
Basic "Live" Observation	8
Capturing a Photo or Video	9-10
Application Menus	10-11
Using the Microscope Tube Off the Base	11
Prepared Slide Specimens	12
Making Your Own Slides	12
Breeding Brine Shrimp	12
Storage	12
Cleaning	12
Warranty	13
FCC Note	13

CONGRATULATIONS ON YOUR PURCHASE !

Thank you for purchasing the Tasco USB Digital Microscope. By combining optics with electronics, this device will help you explore the natural world around you by allowing you to view the microscopic details of plants, animals and minerals on your PC computer monitor.

Before you begin using the USB Microscope, please take the time to read these operating instructions completely, including all special notes and safety guidelines. This will increase your enjoyment of the Microscope, and reduce the risk of personal injury and/or damage to the environment.

Safety Information

This microscope is designed for observing preparations on slides at magnification levels of between 10x to 120x. The enclosed software allows digital recording of the preparations in photo and video format.

Use the USB Microscope only for the purpose described above.

- **This product is intended by the manufacturer for use by persons 13 years of age or older.**
- Keep packing foil and bags out of reach of babies and small children - danger of suffocation!
- The microscope should be connected to computers with USB interface version 1.1 or later. Only use the USB cable supplied. Install the cable in such a way that nobody can trip over it.
- Remove the USB cable from the USB port of the computer when not using the microscope, and prior to any cleaning of the microscope.
- Regularly check all cables and electrical components for any defects or wear. Damaged cables, plugs or other components represent a safety hazard.
- Always position the device on a stable, vibration-free and dry surface.
- Do not expose the device to direct sunlight or heat sources.
- Only use the device inside. Do not use near liquids or in humid rooms.
- Do not use the supplied experimenting materials with objects coming in contact with food. Do not eat, drink or smoke at the place where experiments are carried out.
- **The supplied Eosin dye can cause irritation when coming into contact with the eyes (R36). Do not inhale the dust. If agent comes into contact with the eyes, rinse thoroughly with water and consult a doctor!**
- **Take care when handling the dye, as it can cause discoloration of textiles, furniture and other surfaces.**

Microscope Technical Data

Power supply USB connection version 1.1 or later (5V)

Data transfer USB connection version 1.1 or later

Magnification 10x, 60x,120x selectable

Image resolution RGB 1280 x 1024 pixels

Video resolution RGB 640 x 480 pixels (with connection to USB 2.0 port)

System Requirements (for PC computer)

- Pentium® III 800 MHz or higher is recommended for DV (Digital Video);
- Microsoft® Windows 98 SE, Windows Me, Windows 2000, Windows XP, Windows Vista or Windows 7 (software is not Mac compatible);
- 128 MB RAM (256 MB or higher is recommended)
- 5400 rpm or higher speed hard drive
- CD-ROM or DVD-ROM drive;
- Monitor capable of displaying thousands or millions of colors (1024 x 768 pixels);
- Windows pointing device (mouse, etc.);
- USB 2.0 interface for highest possible video resolution (connection via USB 1.1 allows a maximum resolution of 320 x 240 pixels per frame)

Included Items

- USB Microscope, consisting of the base stand and removable microscope tube
- USB cable
- CD with USB 2.0 PC Camera application/driver (*Windows only*)
- Instruction Manual

(SHOWN BELOW)

- Box of 12 slides (5 prepared and 7 empty slides)
- 4 Sample flasks (gum media (*adhesive*), Eosin dye, sea salt, brine shrimp eggs)
- Storage container and culture dish
- Plastic tweezers

PARTS GUIDE

PARTS GUIDE

- | | |
|--------------------------------------|---------------------------------------------|
| 1. Microscope tube unit | 9. Slide retaining clamps |
| 2. Light button | 10. Microscope stage |
| 3. Ready light | 11. Light: Backlight |
| 4. Control card with image converter | 12. Microscope base (including stand) |
| 5. Snapshot button | 13. Focus knob (on both left & right sides) |
| 6. Magnification dial | 14. USB cable |
| 7. Lens changer | 15. Port for USB cable (rear) |
| 8. Light: Top light | |

ASSEMBLY and SOFTWARE INSTALLATION

Assembly

Remove the Microscope **Tube** (top section, has **Magnification Dial**) and **Base** (lower section, has dual **Focus Knobs**) from the packaging and place the base on a straight, stable and clean surface. Slide the tube from the top into the cradle arms on the base, until the wide section around the **Magnification Dial** rests against the cradle arms (as shown in the parts guide photo). The metal terminals at the rear of the microscope tube should be in contact with the corresponding terminals at the rear of the cradle on the base. These terminals carry power from the **USB Port** to the light in the microscope base.




Software Installation

The microscope's camera will not be recognized by your computer until the driver and application software have been installed on your PC from the included CD, and the computer is restarted. Follow these steps to install the software:

1. Quit any programs that are running on your PC before beginning the installation. Make sure you are logged in to the computer with an administrator account.
2. Place the supplied software CD into the CD or DVD drive. The installation application on the CD will normally open automatically, soon after the CD starts running. *If the CD does not start up automatically, open Windows Explorer and select the drive containing the CD. To start the install menu double click on "AutoRun.exe".*
3. Click "Install PC Cam" on the opening screen (**Fig. 1**). You will get a pop-up window with the message "Please plug in the PC-Cam and click [OK]".
4. Connect the microscope to your PC with the supplied USB cable by connecting the cable first to the microscope (**Fig. 2**) and then to the USB port of your PC. Be sure to connect it to a USB port located directly on the computer, so it will receive power via the USB cable. Do not connect the microscope to an unpowered USB "hub", keyboard USB port, etc. *The microscope should be connected to a computer with ports supporting USB version 1.1 or later. For the highest possible video resolution (640x480), connect to a computer with a USB 2.0 port.*
5. After connecting the microscope, click OK on the pop-up window. The InstallShield Wizard welcome screen will appear (**Fig. 3**). The language for all installation screens is automatically selected based on the language preference setting of your computer. Click on "Next", then Click "Install" on the screen that follows to start the installation (**Fig. 4**). *If you experience a "File Error" popup warning (Fig. 6) at any point during the install process, click "Abort" and eject the CD. You must be logged in to the computer using an administrator account with permission to install software. After logging back in with administrator privileges, reinsert the CD and begin the installation again.*
6. Once the installation has finished, you'll see the "InstallShield Wizard Complete" screen (**Fig. 5**). As it directs, eject the installation disk from the computer's CD drive, then click "Finish" to restart your computer. You have the option of restarting your computer later if you prefer, but you must restart your PC before you begin using the microscope, so it will be properly recognized by the driver.
7. After your computer restarts, you are now ready to begin using the USB Microscope.

Note: If possible, always connect the USB microscope to the same USB port of your PC every time you use it, to avoid additional installation steps and multiple installations of drivers.

SOFTWARE INSTALL SCREENS

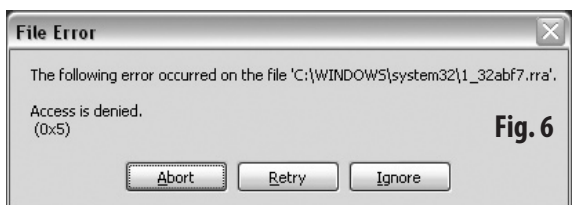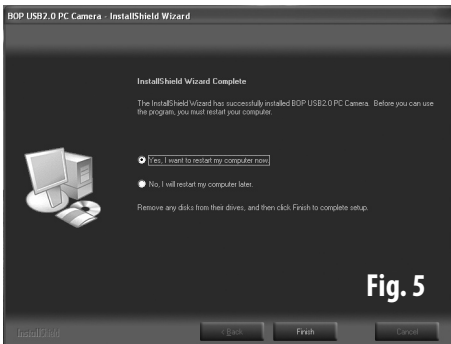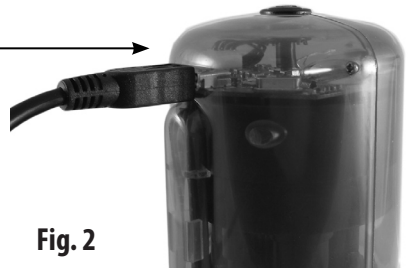USING the MICROSCOPE

Selecting the Lighting

1. Make sure the Microscope is connected to a USB port located directly on the computer, to ensure it will receive adequate power via the USB cable (used for the microscope's lighting, as well as the digital imaging system).
2. Place a slide from the kit under the clips on top of the microscope stage (you can start out with one of the prepared slides).
3. Switch on the microscope's lighting system by pushing the **Light button** one or more times. Your lighting options are:
 - Press 1x: the object is illuminated from above (top lighting). This setting should be used for examining non-transparent objects.
 - Press 2x: the object is illuminated from below (back lighting). This setting should be used for examining transparent objects.
 - Press 3x: the object is illuminated from below and above. Use this setting to show internal and surface structures of semi-transparent objects.
 - Press 4x: Illumination is switched off completely.

Setting the Magnification

The inside of the microscope unit contains 3 optical systems with different magnification factors (10x, 60x, 120x), selected by the **Magnification Dial** at the center of the main tube. The dial can be turned in any direction until it engages. The currently selected magnification is shown on the dial. **Note: Always start your observations on the lowest magnification (10X). This makes it easier to initially center and focus the object. "Finding" the tiny subject on a slide is much more difficult at higher powers, and focusing requires more precision.**

Basic "Live" Observation

1. As described above, make sure the Microscope is connected to your PC via USB, a slide is on the stage ready to view, at least one of the light sources is turned on, and the magnification is set to 10x to start.
2. Launch the application which allows you to view and capture images from the microscope by double clicking the "BOP App" icon on your desktop. *You can also start the application by clicking on the "Camera STI" icon in your Windows system tray (bottom right of screen).*
3. Once the microscope is ready, the **Ready light** comes on. The magnified image of the specimen on the slide now appears in the preview area of the program window.
4. Adjust the image focus by turning the **Focus dial** (on either side).
5. If necessary, move the slide until the subject is fully lit and at the center of the image.
6. You can select a higher magnification if desired. You will normally then need to re-adjust the focus to get a sharp image again.

**BOP App
Icon**

**Camera STI
Icon**

Capturing a Photo or Video

1. In the file menu at the top of the application window, click the File menu and select "Set snapshot file Folder". Choose a folder (for example, "My Pictures") where you would like to save any photos from the microscope you capture with the application. Next, select "Set capture file Folder" from the File menu and choose a folder to store videos you capture with the application.

USING the MICROSCOPE

2. To take a still photo, go to the Capture menu and select "Still Image Capture". This opens a new window showing the captured photo. This window has its own File menu-select Save to store the photo in the folder you choose in the previous step (you can rename the photo, or accept the default file name). Select Exit to close the capture preview window. Note-you can also take a still photo by pressing the Capture button on the front of the microscope tube (just below the green light). However, it may be difficult to avoid handling vibration, resulting in blurry photos-especially above 10x magnification settings.
3. To capture a video clip, select "Start Capture" to begin recording a video clip. Recording will automatically stop when the time limit is up (to change the default time, select "Record time limit" in the Capture menu), or you can manually stop recording video sooner by clicking the "Capture" menu again and selecting "Stop Capture". The .avi video clips are saved to the folder you set up in step #6 above.
4. That's it! When finished using the application, select "Exit" in the File menu. If you want to make adjustments to the still photo or video image quality (this is normally not necessary to obtain good results), explore the options found in the application menus under Option>Video Capture Filter, Option>Video Capture Pin, and Capture>Set Frame rate. These are described in the following Application Menus section.

APPLICATION MENUS:

File

- **Set capture file Folder**-allows you to specify default location for video files (created after selecting "Start Capture" in the Capture menu)
- **Set snapshot file Folder**-allows you to specify default location for still photo files (created after selecting "Still image Capture" in the Capture menu)
- **File index reset**-starts file numbering over with "Still0001" etc. Previous files with same name will be replaced with newer ones.
- **Exit**-quit application

Device: the "USB PC Camera" (your microscope) should be the only device currently using the driver, so normally there is no need to use this menu. However, if you are not getting an image from the microscope (assuming it is connected properly), make sure that "USB PC Camera" is checked in this menu.

- **USB PC Camera**

Option

- **Preview**-leave checked normally. Unchecking and checking again will refresh the "live" image from the camera.
- **Video Capture Filter**-options in the "General Setting" tab allows you to flip the image horizontally ("mirror") or vertically, view the image in black & white, adjust the color saturation, sharpness and brightness and other variables. Note any adjustments you make here will also affect captured still images, not just video clips. The "Special Effect" tab has a variety of entertaining visual effects that can be applied to the microscope image, as well as a digital zoom slider for additional magnification (the "Zoom Enable" box must be checked first). Click "Reset" to restore the original default settings. Click "Apply" to see the Preview image with your current settings. Click OK to keep and switch to your new settings. Click "Cancel" to leave this screen without making any changes.

USING the MICROSCOPE

- **Video Capture Pin**-the video frame rate and output size may be adjusted here if desired to reduce the file size of captured video clips. Normally the frame rate can be left at 30.00 and the output size at the default setting of 1280x1040, unless you find a specific need to change them.

Capture

- **Still image capture**-opens a new window showing the captured still photo, with file menu options allowing you to:
 - Save*- Saves photo to default folder you set in "Set snapshot file Folder". The next default file name is used (to default folder you set in "Set snapshot file Folder"). Files are numbered in a series beginning with "...0001" and continuing unless you use "File index reset" to start over.
 - Save As*- use this instead of "Save" when you want to give your file a name you create (instead of the numbered default name). Using a previous file name again will replace the original, older file with the current captured image.
 - Exit*-select this after you have saved the photo, or use it to cancel if you change your mind or clicked image capture by mistake.
- **Start Capture/Stop Capture**-select "Start Capture" to begin recording a video clip. Recording will automatically stop when the time limit is up (you can set the limit-see below), or you can manually stop recording video sooner by clicking the "Capture" menu again and selecting "Stop Capture".
- **Video format>AVI**-this is already selected, and is the only choice of video format. The captured .avi videos may be viewed with Widows Media Player and most other video software.
- **Record Audio**-this will not affect videos captured by your microscope, so leave it unchecked.
- **Set Frame rate**-Maximum setting is 30 frames per second. Higher frame rates produce smoother video, setting lower rates will generate smaller video files that take up less storage space (but may look "jumpy").
- **Record time limit**-allows you to set the maximum time for each video clip that begins recording when you select "Start Capture". Recording will automatically stop at the end of the time limit (unless you end it early with "Stop Capture").
- **Enable BPC> Y or UV**-This should be left on the default settings, with both checked.

Help

- **About Cam App**-displays software version number

Using the Microscope Tube "Off the Base" (Optional)

The **Microscope main tube** (top section) may be used independently of the base. This provides the option of examining details on larger subjects or objects that cannot physically fit on the slide stage area of the base. To use the tube away from the base, simply reverse the assembly process and carefully slide the tube straight up away from the base. **Make sure the USB cable remains attached, and that you have enough slack in the cable to safely move the tube towards your subject without risk of damage to the computer.** As the focus knobs are part of the base, you will need to focus by slowly moving the tube closer to or away from the subject until it appears sharp (using the 10X magnification setting is recommended). **Be careful not to touch the subject with the surface of the lens to avoid scratches or other damage.** Note that only the top light is available when using the tube alone.

USING the ACCESSORIES

Your Tasco USB Digital Microscope is a versatile instrument. To start off beginning science enthusiasts, the set contains also some accessory materials to help them perform basic observation without the need for additional specialized equipment.

Prepared Slide Specimens

The following ready-made slides are part of the contents of the box and can be analyzed under the microscope without further preparation:

- Apple
- Plant pollen
- Leaf of water thyme (*Hydrilla verticillata*)
- Leaf of broad bean (*Vicia faba*)
- Onion skin (*Allium cepa*)

Making Your Own Slides

The following accessories have been provided for preparing your own slides:

- **Slide:** Using a sharp blade cut a very fine sample from the material to be examined and place it on the clear slide.
- **Adhesive** (Rubber Arabicum): This substance is water-soluble and is used to attach preparations to the slide.
- **Eosin coloring:** Eosin is a strong dye for coloring preparations, making it easier to see details (such as cell walls or cell nuclei). The use of minute amounts together with a drop of water suffices.

Caution!

The supplied Eosin dye can cause irritation when coming into contact with the eyes. Do not inhale the dust! If agent comes into contact with the eyes, rinse thoroughly with water and consult a doctor! Use care when handling the dye. It can cause discoloration of textiles, furniture and other surfaces.

Breeding Brine Shrimp

An interesting experiment is observing the development of brine shrimp (*Artemia Salina*) under the microscope. Multiple sources of information describing in detail the process of hatching and raising brine shrimp are readily available on the internet.

To observe the animals always place them in a salt-water droplet on the slide or place them under the microscope in a Petri dish filled with salt water.

Storage

When not in use, store your USB Digital Microscope in a dry location free from dust. If the device is not used for some time it is recommended that it be returned to the original packaging for secure storage and also transportation.

Cleaning

Your USB Microscope is a high-quality optical device. Avoid contact with dust and humidity. Avoid fingerprints on all optical surfaces. First remove dirt or dust using a soft brush. Next, clean the soiled area with a soft, lint-free cloth (e.g. linen) and some alcohol. Do not use any other solvents or aggressive or abrasive cleaning agents as these could damage the surface.

ONE YEAR LIMITED WARRANTY

Your Tasco® product is warranted to be free of defects in materials and workmanship for one year after the date of purchase. In the event of a defect under this warranty, we will, at our option, repair or replace the product, provided that you return the product postage prepaid. This warranty does not cover damages caused by misuse, improper handling, installation, or maintenance provided by someone other than a Tasco Authorized Service Department.

Any return made under this warranty must be accompanied by the items listed below:

- 1) A check/money order in the amount of \$10.00 to cover the cost of postage and handling
- 2) Name and address for product return
- 3) An explanation of the defect
- 4) Proof of Date Purchased
- 5) Product should be well packed in a sturdy outside shipping carton, to prevent damage in transit, with return postage prepaid to the address listed below:

IN U.S.A. Send To:

Tasco Outdoor Products
Attn.: Repairs
9200 Cody
Overland Park, Kansas 66214

IN CANADA Send To:

Tasco Outdoor Products
Attn.: Repairs
25A East Pearce Street, Unit 1
Richmond Hill, Ontario L4B 2M9

For products purchased outside the United States or Canada please contact your local dealer for applicable warranty information. In Europe you may also contact Tasco at:

Tasco Germany GmbH
European Service Centre
Mathias-Brüggen-Str. 80
D-50827 Köln
GERMANY
Tel: +49 221 995568-0
Fax: +49 221 995568-20

This warranty gives you specific legal rights.

You may have other rights which vary from country to country.

©2010 B.O.P.

FCC NOTE

This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

Shielded interface cable must be used with the equipment in order to comply with the limits for a digital device pursuant to Subpart B of Part 15 of FCC Rules.

Specifications and designs are subject to change without any notice or obligation on the part of the manufacturer.

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Disposal of Electric and Electronic Equipment

(Applicable in the EU and other European countries with separate collection systems)

This equipment contains electric and/or electronic parts and must therefore not be disposed of as normal household waste. Instead, it should be disposed at the respective collection points for recycling provided by the communities. For you, this is free of charge.

If the equipment contains exchangeable (rechargeable) batteries, these too must be removed before and, if necessary, in turn be disposed of according to the relevant regulations (see also the respective comments in this unit's instructions).

Further information about the subject is available at your community administration, your local waste collection company, or in the store where you purchased this equipment.

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